

SHIMSEN Mini-Filter Vial




SHIMSEN Mini-Filter Vial simplify your workflow and reduce waste generated in the lab by replacing four different components with one Mini-Filter Vial.

- Replaces syringe, syringe filter, vial, and cap;
- Waste and cost reduction;
- Includes visual indication that the sample has been filtered;
- Minimizes instrument downtime due to unfiltered samples;





■ How to use?

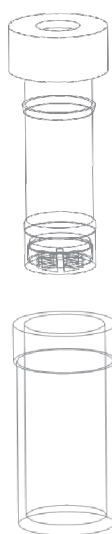


Fill

Use a pipette to add sample to the fill line.

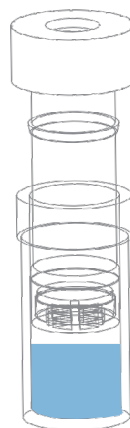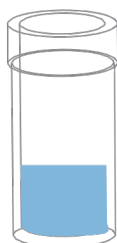

Cover

Twist gently to ensure a secure seal.

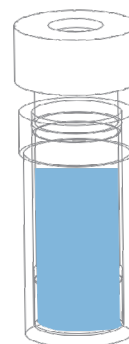

Plunge

To filter the sample, press the plunger slowly.

SHIMSEN Mini-Filter Vial

P01-P03



380-06211-01



380-06211-02



380-06211-03



380-06211-04

Ref.No.	380-06211-01	380-06211-02	380-06211-03	380-06211-04
Description	Nylon 0.22 µm Clear Filter Vial,pp	Nylon 0.45 µm Clear Filter Vial,pp	PTFE 0.22 µm Clear Filter Vial,pp	PTFE 0.45 µm Clear Filter Vial,pp
100 pcs/pk.				



380-06211-05



380-06211-06



380-06211-07



380-06211-08

Ref.No.	380-06211-05	380-06211-06	380-06211-07	380-06211-08
Description	PES 0.22 µm Clear Filter Vial,pp	PES 0.45 µm Clear Filter Vial,pp	PVDF 0.22 µm Clear Filter Vial,pp	PVDF 0.45 µm Clear Filter Vial,pp
100 pcs/pk.				



380-06211-09



380-06211-10



380-06211-11



380-06211-12

Ref.No.	380-06211-09	380-06211-10	380-06211-11	380-06211-12
Description	Nylon 0.22 µm Amber Filter Vial,pp	Nylon 0.45 µm Amber Filter Vial,pp	PTFE 0.22 µm Amber Filter Vial,pp	PTFE 0.45 µm Amber Filter Vial,pp
100 pcs/pk.				

SHIMSEN Mini-Filter Vial

SHIMSEN Mini-Filter Vial

P01-P03



380-06211-13



380-06211-14



380-06211-15



380-06211-16

Ref.No.	380-06211-13	380-06211-14	380-06211-15	380-06211-16
Description	PES 0.22 μm Amber Filter Vial,pp	PES 0.45 μm Amber Filter Vial,pp	PVDF 0.22 μm Amber Filter Vial,pp	PVDF 0.45 μm Amber Filter Vial,pp
100 pcs/pk.				



380-06211-17



380-06211-18



380-06211-19



380-06211-20

Ref.No.	380-06211-17	380-06211-18	380-06211-19	380-06211-20
Description	Ultra-clean PTFE 0.22 μm Clear Filter Vial,pp	Ultra-clean PTFE 0.45 μm Clear Filter Vial,pp	Ultra-clean PTFE 0.22 μm Amber Filter Vial,pp	Ultra-clean PTFE 0.45 μm Amber Filter Vial,pp
100 pcs/pk.				

Product size : Diameter 12 mm, Height : 34mm

Sterile: No

Filtration Area: 0.28cm²

Filling vol/Dead vol : 480ul/30ul

Max Pressure: 1.5 bar (21.6 psi)

Max operation temp: 121 °F (50 °C)

Standard particle retention efficiency ≥ 95%



Shimadzu (Shanghai) Global Laboratory Consumables Co.,Ltd.

www.sglc.shimadzu.com.cn

www.shimadzumall.com

Contact: contact@sglc.shimadzu.com.cn

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. See <http://www.shimadzu.com/about/trademarks/index.html> for details.

Third party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they are used with trademark symbol "TM" or "®".

Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.