


# Small Molecule Pharmaceutical Supplies Selection Guide

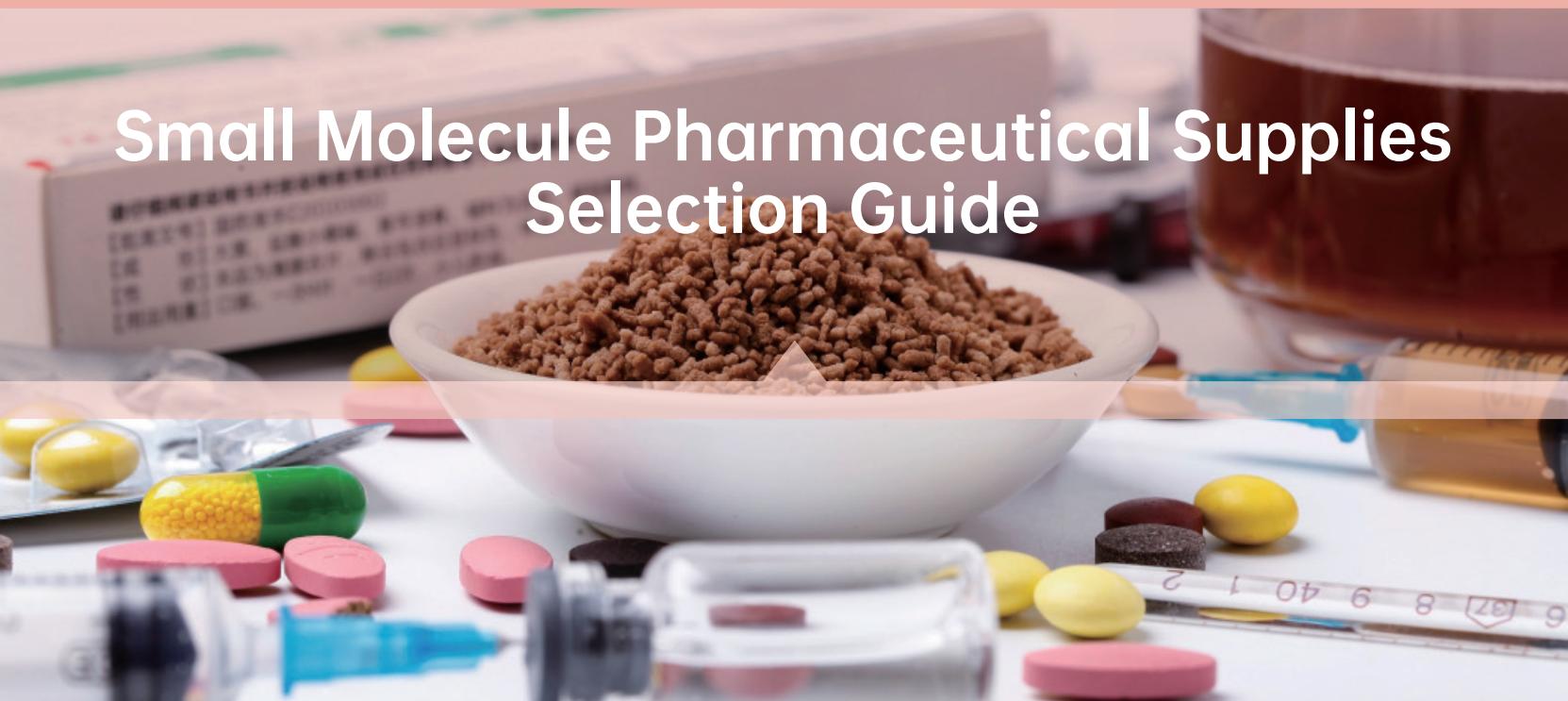

# SGLC offers "one-stop" consumables solutions

## You are not alone with SGLC

### “Preface

SGLC is providing this Small Molecule Pharmaceutical Supplies Selection Guide based on the needs of Pharmaceutical Analysis.



# CONTENTS



1 Related Substances

2 Genotoxic Impurities

3 ASSAY

4 Residual Solvents

5 Dissolution Test

6 Excipients&Packing Materials

7 Preparative Scale LC Column


8 Featured Solutions

# Related Substances

Separation is the primary concern for related substances analysis.

## Shim-pack Scepter Series




- The next generation organic silica hybrid based columns!
- A wide range of LC conditions!
- Best choice of Related substances analysis

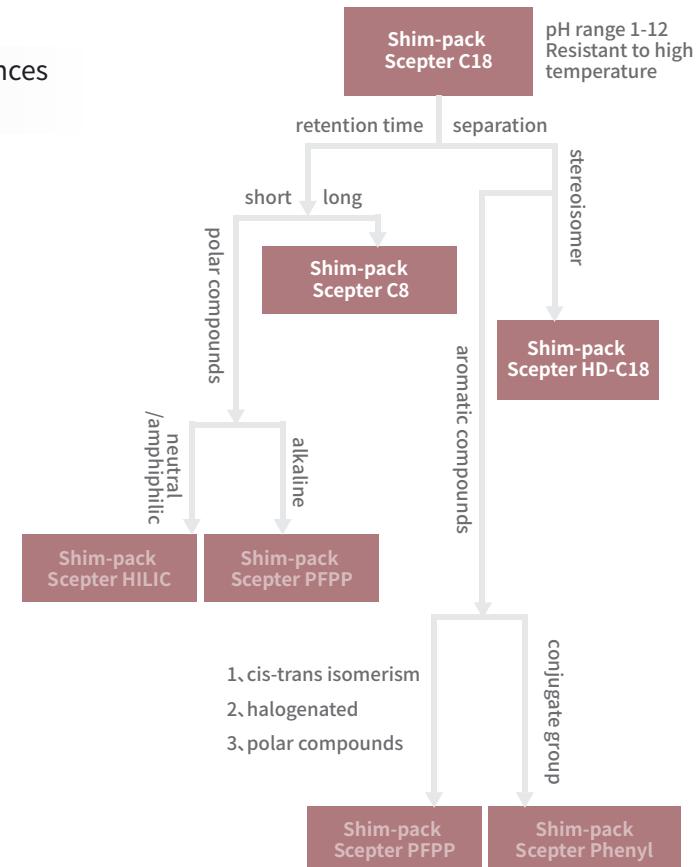


# Reversed phase


Classification	Drug	LC Column	Feature	Classification	Drug	LC Column	Feature
Antibiotics	Azithromycin, Penicillin Sodium, Cefaclor, Cefazolin Sodium, Cefoperazone Sodium	Shim-pack Scepter C18 (4.6mm×250mm, 5μm, PN: 227-31020-06)	Good separation and reproducibility.	Anti-inflammatory	Aceclofenac	Shim-pack Scepter C18 (4.6mm×250mm, 5μm, PN: 227-31020-06)	Perfect separation and peak symmetry.
	Sulbactam Sodium, Flucloxacillin Sodium	Shim-pack Scepter C18 (4.6mm×150mm, 3μm, PN: 227-31016-05)	Good separation of impurities and API.		Dexibuprofen	Shim-pack Scepter C18 (4.6mm×150mm, 3μm, PN: 227-31016-05)	Baseline separation of isomeric impurities.
Hormone	Rednisone Acetate, Hydrocortisone, Cortisone Acetate, Hydrocortisone Acetate, Ethinyl Estradiol, Triamcinolone Acetate	Shim-pack Scepter C18 (2.1mm×100mm, 1.9μm, PN: 227-31012-05)	UHPLC method with good separations& fast.	SHIMADZU Artemisinin-based RIPPLE PACKED C18	Chlortetracycline	Shim-pack Scepter C18 (4.6×100 mm, 3 μm, PN: 380-31016-04)	Symmetrical peaks & good separation.
	Dexamethasone Acetate	Shim-pack Scepter C18 (4.6mm×250mm, 5μm, PN: 227-31020-06)	Baseline separation & good peak shapes.		Dihydroartemisinin, Artesunate	Shim-pack Scepter C18 (4.6×100 mm, 3 μm, PN: 380-31016-04)	Good separation & perfect peak shapes.
Antihypertensive	Rbesartan, Captopril	Shim-pack Scepter C18 (4.6mm×150mm, 3μm, PN: 227-31016-05)	Good separation& perfect peak shapes.	Antivirals	Adefovir	Shim-pack Scepter C18 (4.6mm×150mm, 5μm, PN: 227-31020-05)	Baseline separation.
	Nifedipine	Shim-pack Scepter C18 (2.1mm×100mm, 1.9μm, PN: 227-31012-05)	Separation Rs>1.5.		Valaciclovir Hydrochloride	Shim-pack Scepter C18 (2.1mm×100mm, 1.9μm, PN: 227-31012-05)	UHPLC method & baseline separation.
	Nimodipine	Shim-pack Scepter HD-C18 (4.6mm×250mm, 5 μm, PN: 227-31024-06)	Baseline separation & perfect peak shapes.	Antihyper-glycemic	Gliquinone	Shim-pack Scepter C18 (2.1mm×100mm, 1.9μm, PN: 227-31012-05)	UHPLC method.
	Perindopril Indapamide	Shim-pack Scepter C8 (4.6mm × 250mm, 5μm, PN: 227-31041-06)	Good separation & symmetric peaks.	Hyperuricemia/gout	Febuxostat	Shim-pack Scepter HD-C18 (4.6mm×250mm, 5 μm, PN: 227-31024-06)	The baseline separation; good peak shapes and reproducibility.
Anti-tumor	Flutamide	Shim-pack Scepter C18 (4.6mm×150mm, 3μm, PN: 227-31016-05)	Perfect peak shape & good separation.	Muscle relaxant	Baclofen	Shim-pack Scepter C18 (4.6 mm×250 mm, 5μm, PN: 227-31020-06) Shim-pack Scepter C18 (2.1 mm×100 mm, 1.9μm, PN: 227-31012-05)	Baseline separation, good retention of the main peak.
	Letrozole	Shim-pack Scepter C18 (4.6mm×150mm, 3μm, PN: 227-31016-05)	UHPLC method achieves good baseline separation of the 4.4min main peak from isomeric.	Antimicrobial	Terbinafine Hydrochloride	Shim-pack Scepter C18 (2.1mm×100mm, 1.9μm, PN: 227-31012-05)	Good separation & perfect peak shapes.

# Related Substances

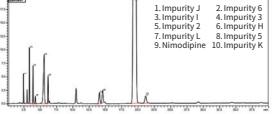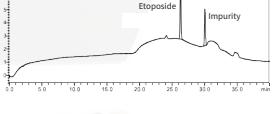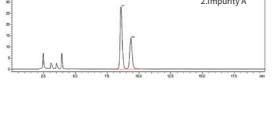Drug	LC column	Feature	Chromatogram
Nimodipine	Shim-pack Scepter Phenyl-120 (4.6 mm × 250 mm, 5 µm, PN:227-31071-06)	Good peak shapes; separation meets the requirements.	
Etoposide	Shim-pack Scepter Phenyl-120 (4.6 mm × 250 mm, 5 µm, PN:227-31071-06)	SHIMADZU Separation and retention time were in accordance with the requirements of pharmacopoeia.	
Levocanдин	Shim-pack Scepter Phenyl-120 (4.6 mm × 250 mm, 5 µm, PN:227-31071-06)	Symmetric peaks & good separation.	

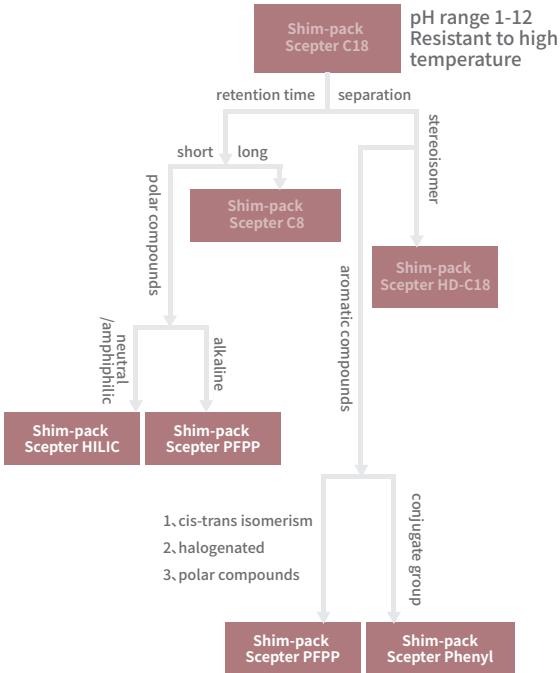

# "Magic tool" for Related Substances Analysis-Ghost Trap DS

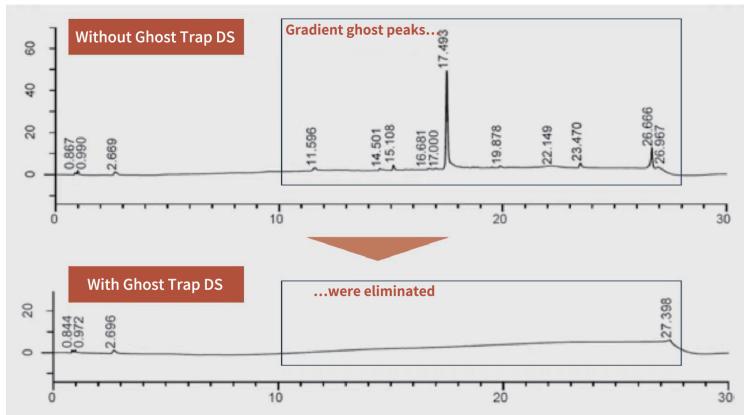

Column : ODS column  
 Mobile Phase : A) 25 mmol/L Phosphate (Potassium) buffer solution (pH 4.0) /Acetonitrile = 9/1  
 B) Water/Acetonitrile = 1/9  
 Flow Rate : 0.65 mL/min  
 Col. Temp : 45 °C  
 Detection : UV 210 nm

Ghost Trap DS\* was co-developed with Daiichi Sankyo Co., Ltd.  
 Higher trapping efficiency and longer life!


\* DS: Abbreviation of Daiichi Sankyo (D) and Shimadzu (S)



Ghost Trap DS

Economic type

TIPS.

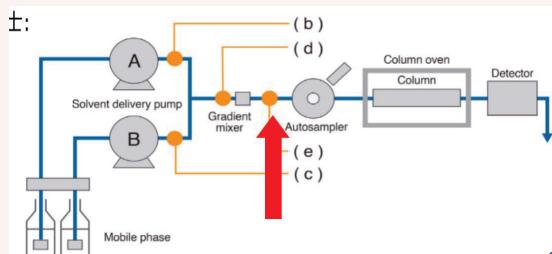

- Installation position: (e). Between the injector and mixer. Only capture the ghost peaks from the mobile phase, pipeline, system; no effect on the sample.
- Not apply to ion-pair reagent.




# Genotoxic Impurities

Better separation, higher sensitivity!

USP <1469> specifies detection methods for nitrosamine-based toxic impurities that include HPLC-MS/MS and GC-MS/MS.

## Application 1 Quantification of Nitrosamines (LCMS/MS)

Shim-pack Velox C18 -- High sensitivity & fast

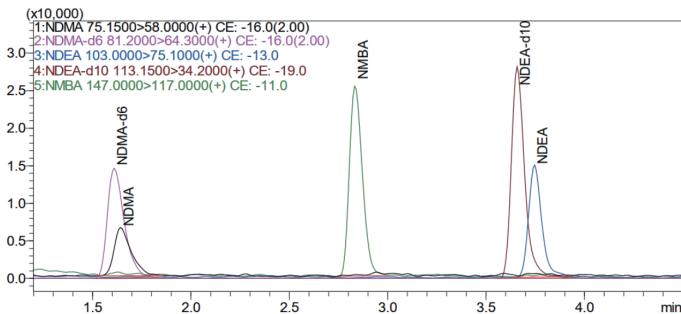

### LC conditions :

Column: Shim-pack Velox C18 (2.1mm × 100 mm, 1.8µm, PN: R227-32007-03)

Eluent: A-0.1% formic acid; B-methanol

Needle wash mode: before and after injection

Flow rate: 0.3 mL/min


Temperature: 40 °C

Injection volume: 20 µL

Elution mode: gradient elution

## Application 2 Structural characterization of Nitrosamines (LCMS-QTOF)

Shim-pack Scepter C18 -- Excellent separation of impurities near the main peak

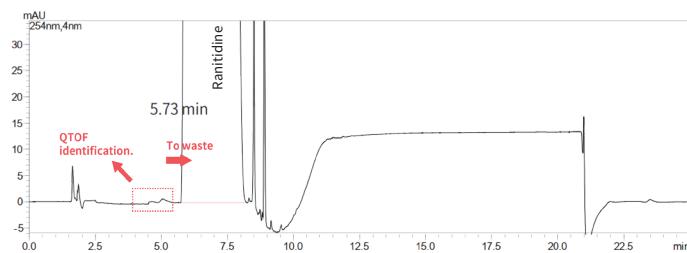

### LC conditions :

Column: Shim-pack Scepter C18 (4.6mm × 150mm, 5.0µm, PN: 227-30017-07)

Eluent: A-0.1% formic acid; B-methanol

Flow rate: 1.0 mL/min


Injection volume: 5 µL

Temperature: 40 °C

Elution mode: gradient elution

### Application 3 Quantification of Nitrosamines (HS-GC-MS)

**SH-I-624Sil MS** -- Excellent linearity and signal-to-noise ratio!

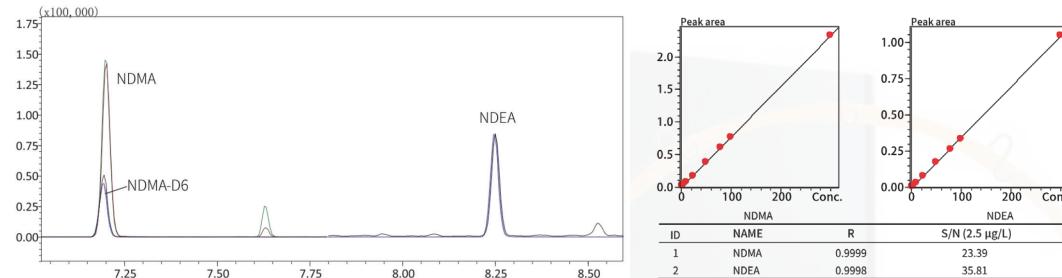

#### GC conditions :

Column: SH-I-624Sil MS (60 m × 0.32 mm × 1.8 µm PN:221-75963-60)

Temperature program: 40°C(1 min)\_40°C/min\_180°C\_20°C/min\_300°C(2 min)

Injection volume: 2 µL

Carrier gas control mode: constant linear velocity (36.1 cm/sec) Inlet mode: non-split injection Ion source temperature: 230°

Interface temperature: 300°C Detector voltage: tuned voltage +0.6kV Acquisition method: MRM

### Application 4 Quantification of Methanesulfonates (HS-GC-MS)

**SH-PolarWax** -- Bonded Crosslink™ polyethylene glycol stationary phase, low loss ensures column lifetime

#### Headspace conditions

Equilibration temperature: 60°C; Equilibration time: 30 min; Sample pathway temperature: 100°C.

#### GCMS conditions

Column: SH-PolarWax Column (30m × 0.25mm × 0.25 µm, PN:227-36305-02)

Temperature program: 40°C(1 min)\_10°C/min\_130°C\_40°C/min\_230°C(8 min)  
Carrier Gas: He, constant flow mode.


Split ratio: 20:1

Ionization mode: EI

Ion source temperature: 220°C

Interface temperature: 220°C

Flow rate: 0.8 mL/min to 2 mL/min after 10 min.

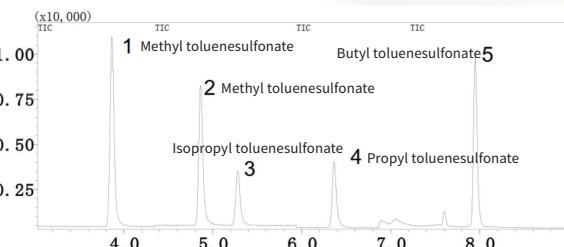

# ASSAY

## ASSAY

### ■ ShimNex UP Series

- Silica-based, multi-specifications!
- High column-to-column performance reproducibility!
- Best choice for ASSAY!




		ShimNex UP Series						
Ligand Type	Particle Size	Pore Size	LD.:mm					
			2.1	3.0	4.0	4.6	10.0	21.2
C18	1.8μm	100 Å	●	●	○	○	○	○
	2.5μm		●	●	○	●	○	○
	3.5μm		●	●	●	●	○	○
	5μm		●	●	●	●	●	●
	10μm		○	○	●	●	●	●
C8	1.8μm	100 Å	●	●	○	○	○	○
	2.5μm		●	●	○	●	○	○
	3.5μm		●	●	●	●	○	○
	5μm		●	●	●	●	●	●
	10μm		○	○	●	●	●	●
C4	1.8μm	100 Å	●	●	○	○	○	○
	2.5μm		●	●	○	●	○	○
	3.5μm		●	●	●	●	○	○
	5μm		●	●	●	●	●	●
	10μm		○	○	●	●	●	●

Length : 50mm, 100mm, 150mm, 250mm

注： ● :available; ○ :unavailable

### Vitamin D2 & D3

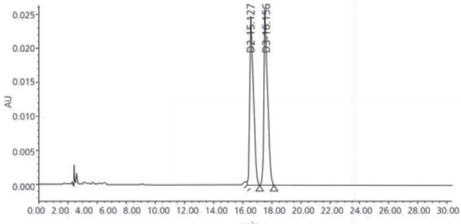

#### LC conditions:

**Column:** ShimNex UP C18, 4.6x250mm, 5 $\mu$ m (PN:380-01231-49)

**Eluent:** methanol:water=98:2


**Flow rate:** 1.0 mL/min

**Temperature:** 40°C

**Detection:** UV 265nm

**Injection volume:** 10  $\mu$ L

### Vitamin E & D

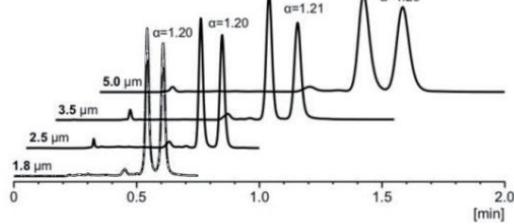

#### LC conditions:

**Column:** ShimNex UP C4, 2.1x50mm, 1.8 $\mu$ m or 4.6x50mm,  $\geq$ 2.5 $\mu$ m


**Eluent:** acetonitrile

**Flow rate:** 5.0  $\mu$ m: 1.0 mL/min; 3.5  $\mu$ m: 1.5 mL/min; 2.5  $\mu$ m: 2.0 mL/min; 1.8  $\mu$ m: 0.6mL/min

**Temperature:** 20°C

**Detection:** UV 215nm

### Amino acids

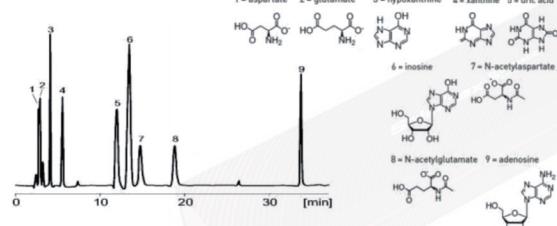

#### LC conditions:

**Column:** ShimNex UP C18, 4.6x250mm, 5 $\mu$ m (PN:380-01231-49)


**Eluent:** 28 mM tetrabutylammonium hydroxide, 25 mM potassium dihydrogen phosphate, 1.25% methanol (pH 7)

**Flow rate:** 1.0 mL/min

**Temperature:** 23°C

**Detection:** UV 210nm

### Norfloxacin

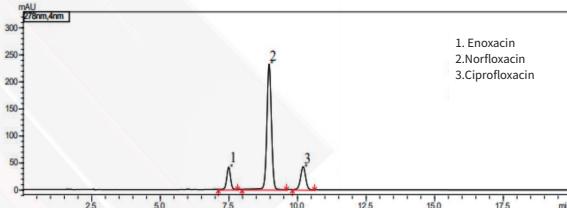

#### LC conditions:

**Column:** ShimNex UP C18, 4.6×250 mm, 5  $\mu$ m (PN:380-01231-49)

**Eluent:** 0.025 mol/L Phosphoric acid (adjusted to pH 3.0 with triethylamine)-acetonitrile=87:13

**Flow rate:** 1.5 mL/min

**Temperature:** 40°C

**Detection:** UV 278nm

**Injection volume:** 20  $\mu$ L

# Analysis of Residual Solvents

## Residual Solvents

### SH-I-624Sil MS


- Mid-polarity Crossbond™ silarylene phase.
- Ultra-inert & high-thermal stability!
- Best choice for residual solvent analysis!



### List of GC Columns for Residual Solvent Analysis

Category	Columns	S/N No.
Non-polar (-1)	SH-I-1MS Cap. Column, 30mx0.32mmx1μm	R227-36010-01
Medium polarity (-624)	SH-I-624Sil MS Cap. Column, 30mx0.32mmx1.8μm	R227-36077-01
Polar (WAX)	SH-PolarWax Column, 30mx0.25mmx0.25μm SH-PolarWax-MS, 30mx0.25mmx0.25μm	R227-36305-02 R227-36322-01
Columns for amine analysis	SH-Volatile Amin Cap. Column, 30mx0.32mmx5μm	R227-36326-01

### Analysis of Class 1 & Class 2 Residual Solvents

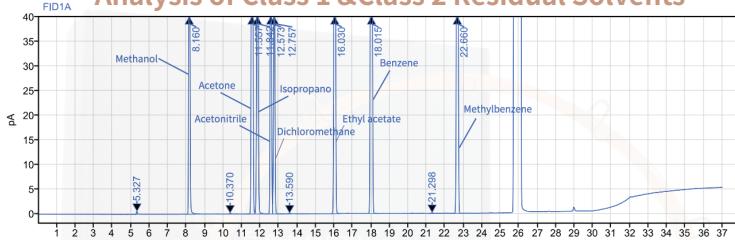

Column: SH-I-624Sil MS, 30 m × 0.53 mm × 3.0 μm (PN: R227-36078-01)

Injection method: Headspace injection

Split ratio: 5: 1

Equilibrium temperature: 80 °C

Inlet temperature: 250 °C


Quantification loop temperature: 100 °C

GC cycle time: 45 min

GC cycle time: 45 min

Temperature program: 35 °C (hold 5 min) to 160 °C at 5 °C/min (hold 5 min) to 240 °C at 40 °C/min (hold 5 min)

### Analysis of Amines

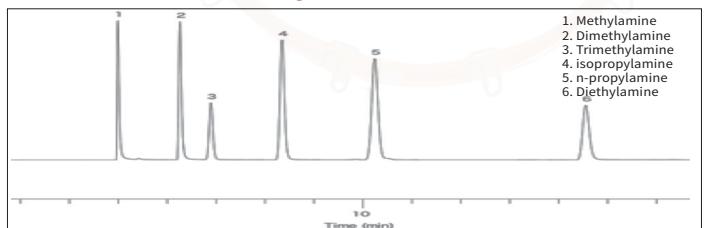

System: GC-FID

Column: SH-Volatile Amin (60 m × 0.32 mm × 5.00 μm)

Temperature: 60°C

Carrier gas: He 150kPa



Split ratio: 1:50

# Dissolution Test

## ■ Shim-pack Scepter Series

- Ultra-high tolerance, fast, easy to handle complex sample matrices!
- The best choice for analyzing dissolved samples!

## In Vitro Dissolution Study of Ambroxol Hydrochloride Tablets in Four Dissolution Medias

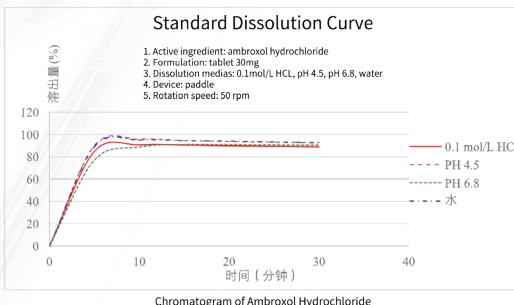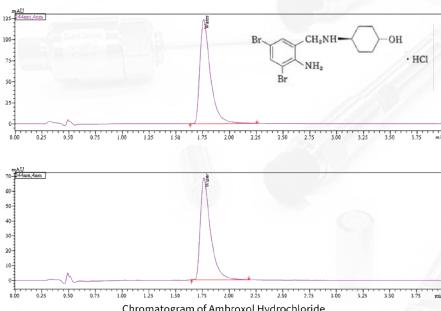

### LC conditions:

**Column:** Shim-pack Scepter C18 (4.6mm×50mm, 5μm, PN: 227-31020-02)

**Eluent:** A- 0.01mol/L Ammonium dihydrogen phosphate B: acetonitrile, A/B = 1:1(v/v)

**Flow rate:** 0.4 mL/min

**Temperature:** 40°C

**Detection:** 244nm

**Injection volume:** 5μL

The conditions of dissolution determination were operated according to the draft of generic drug consistency evaluation of Amoxicillin Hydrochloride Tablets drafted by National Medical Products Administration.



## TIPS.



Reasons for unstable retention time: column contamination, insufficient column equilibration time, selective volatilization due to different nature of mobile phase, unstable column temperature



# Pharmaceutical Excipients & Packing Materials

## ■ SH Series GC Columns

- Excellent chromatographic performance, reproducibility & lifetime!

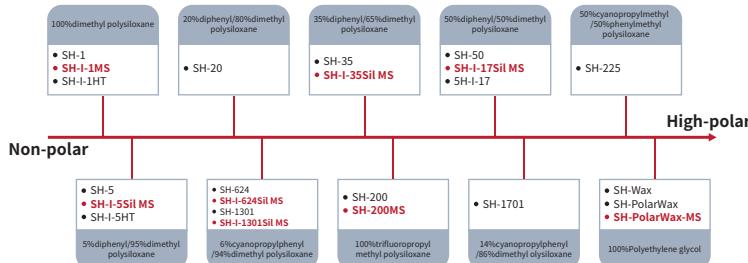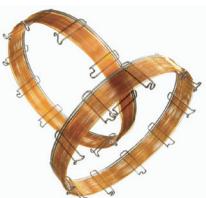

Cases	GC Column	Features	Chromatograms
<b>Determination of ethylene oxide (EO) in pharmaceutical packaging materials (pre-filled syringes)</b>	SH-624 30 m×0.25 mm × 1.4 µm PN:R221-75863-30; SH-I-624Sil MS 30 m×0.32 mm × 1.8 µm PN:R227-36077-01	Ultra-high inertness and low loss, temperature resistant to 300/320°C, unique selectivity for volatile organic compounds with boiling points below 200°C, such as ethylene oxide, propylene oxide, epichlorohydrin, etc. EO peak is baseline separated from the adjacent impurity acetaldehyde peak in the spectrum.	
<b>Glycol, diethylene glycol, triethylene glycol in Polysorbate 80</b>	SH-PolarWax 30 m×0.25 mm × 0.25 µm PN:227-36305-02	Less interference between ethylene glycol and propylene glycol; better peak shapes!	
<b>Fatty acid composition of lipid substances for pharmaceutical and excipients</b>	SH-Wax 30 m×0.53 mm × 1 µm PN:221-75899-30	Determination of 36 fatty acids, meets the system suitability requirements of the fatty acid composition determination of the regulation.	

# Preparative Scale LC Column

## Shim-pack Scepter PREP series

- A wide range of LC conditions & multiple phases!
- High column-to-column performance reproducibility
- Best choice of Preparative columns




**Flow rate equation (Same with Loading) :**


$$\text{Prep flow rate} = \text{Analytical flow rate} \times \frac{(\text{Prep I.D.})^2}{(\text{Analytical I.D.})^2}$$

Lab Scale → Production Scale										
	Loading (mg)	Lab Scale	10-50	50-100	100-1000	1g-10g	10g -100g	≥100g		
High Efficiency, Cost, Purserse ↓ Low	Particle Size (μm)	I.D.(mm)	4.6	10	20	30	50	100	200	≥300
	5	●	●	●	●	●	●	●	●	●
	10	●	●	●	●	●	●	●	●	●
	10-20	●	●	●	●	●	●	●	●	●
	15-30	●	●	●	●	●	●	●	●	●
	50				●	●	●	●	●	●


● Most appropriate    ● Appropriate    ● Depending on purpose

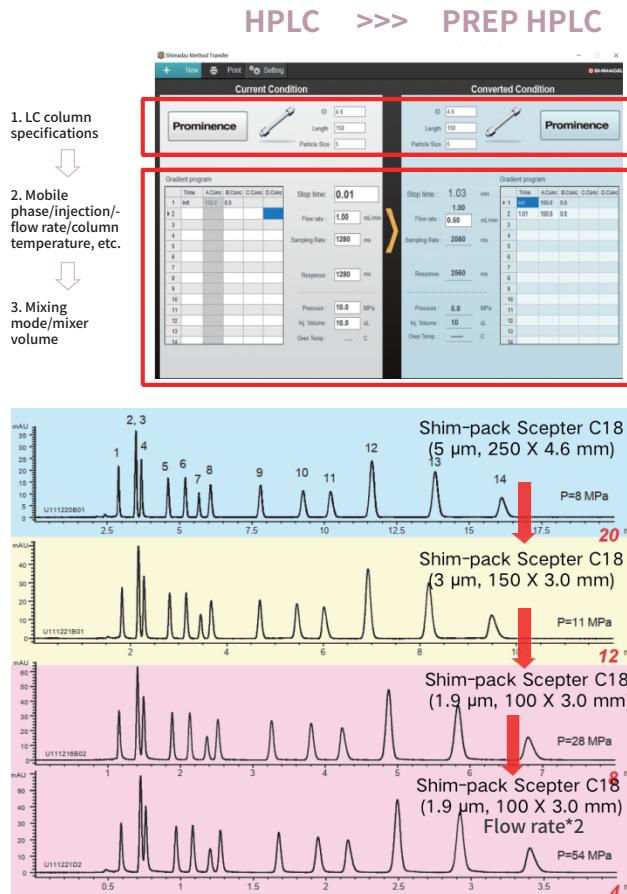

# Featured Solution- desalted by 2D chromatography & identification of cephalosporin polymers

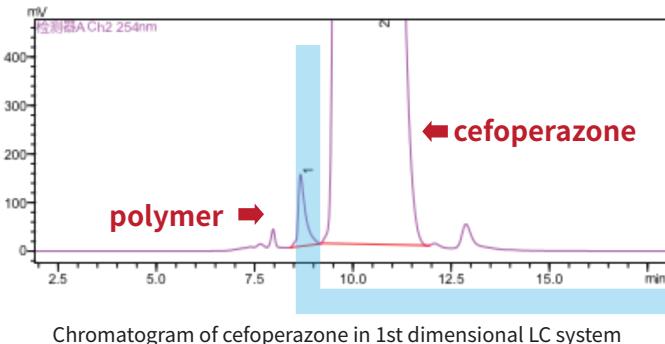

Chromatogram of cefoperazone in 1st dimensional LC system

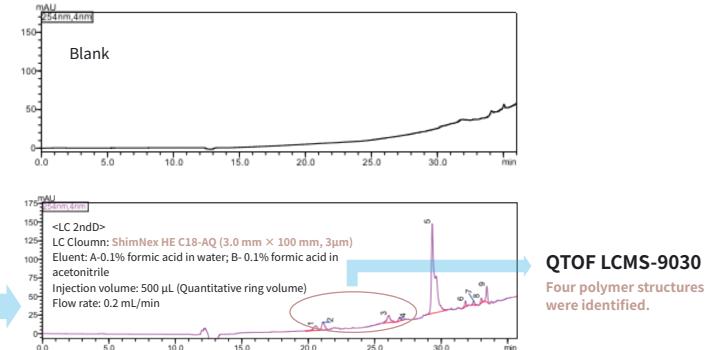

TIPS.

## SHIMSEN Ankylo-SEC

Fast & high column efficiency;  
excellent batch reproducibility.



## ShimNex HE C18-AQ

Homogeneous particle size, rapid  
elution under the premise of  
ensuring the separation.

- ▶ Non-volatile mobile phase additives are not allowed to enter the mass spectrometry detector.


# Featured Solution- column temperature adjustment

Improvement of separation and peak shapes by increasing column temperature

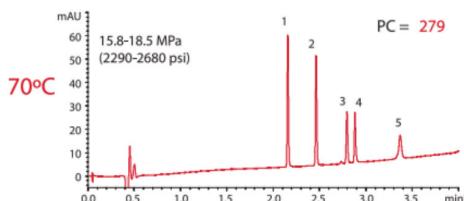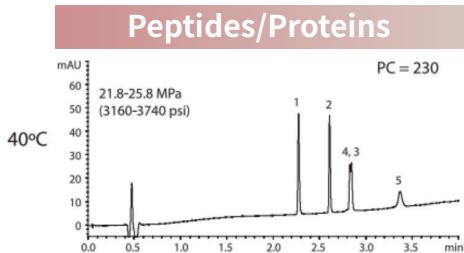


Analytes	MW	Peak width <sub>0.5h</sub> (min)	
		40°C	70°C
<b>Condition A</b>			
1. Oxytocin	1,007	0.017	0.014
2. Leu-Enkephalin	556	0.015	0.015
3. $\beta$ -Endorphin	3,465	-	0.016
4. Insulin	5,733	-	0.015
5. $\beta$ -Lactoglobulin A	18,400	0.043	0.030

### Macrolide Antibiotic

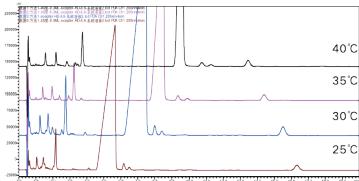

### Ribonucleic acid

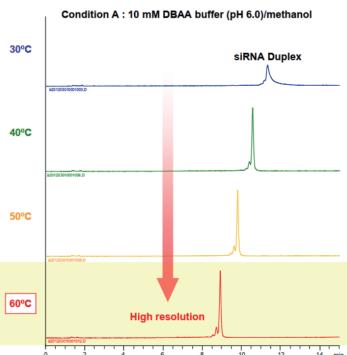


### Shim-pack Scepter C18

Organic silica hybrid based column  
pH range 1-12  
High temperature resistance



Stability at high pH  
(pH 11.5, 40°C)

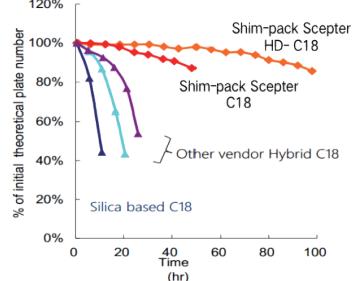

TIPS.

Be sure to use the column within the specified conditions.

# 化药相关消耗品推荐

**Vials**



SHIMSEN Vial  
Laboratory General Vials



SHIMSEN ARC Vial  
Economic choice



SHIMSEN Low adsorption  
Low adsorption PP material



Labtotal Vial  
LC/MS and GC/MS certification report  
Cover gasket pre-installation

**Needle Filter**



HPTFE  
For aquatic and organic



PTFE  
For organic



PES  
For Aquatic

**Other Consumables**



Safety caps



Solvent bottle



Pipette

## Equipment



Nitrogen generator for LC-MS



Air compressor for GC



Hydrogen generator for GC



Air compressor for AA



Cooling circulating water unit for AA

# Book Online

岛津消耗品全套方案 一站式解决!

## 行业解决方案



## 产品选型卡





**SHIMADZU**

**Shimadzu (Shanghai) Global Laboratory  
Consumables Co.,Ltd.**

www.sglc.shimadzu.com.cn  
www.shimadzumall.com  
Contact: contact@sglc.shimadzu.com.cn

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu.  
See <http://www.shimadzu.com/about/trademarks/index.html> for details.

Third party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they are used with trademark symbol "TM" or "®".

Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.